# Levenhuk MED 5M Microscope Digital Camera with 9.4" LCD Screen

- EN User manual
- BG Ръководство за потребителя
- CZ Návod k použití
- DE Bedienungsanleitung
- ES Guía del usuario
- HU Használati útmutató
- IT Guida all'utilizzo
- PL Instrukcja obsługi
- PT Manual do usuário
- RU Инструкция по эксплуатации

*Наслади се отблизо  
Radost zaostřit  
Zoom ran und hab Fun!  
Amplíe y disfrute  
Kellemes nagyítást!  
Ingrandisci il divertimento  
Radość przybliżania  
Dê um zoom na sua emoção  
Приближает с удовольствием*

**levenhuk**  
Zoom&Joy

<b>EN</b>	Levenhuk MED 5M Microscope Digital Camera with 9.4" LCD Screen	<b>3</b>
<b>BG</b>	Цифрова камера за микроскоп Levenhuk MED 5M с 9,4" течнокристален екран	<b>14</b>
<b>CZ</b>	Levenhuk MED 5M Digitální Fotoaparát Mikroskopu s 9,4" LCD obrazovkou	<b>25</b>
<b>DE</b>	Levenhuk-Mikroskopdigitalkamera MED 5M mit 9,4-Zoll-LCD-Bildschirm	<b>36</b>
<b>ES</b>	Cámara digital de 5M para microscopio Levenhuk MED con pantalla LCD de 9,4"	<b>47</b>
<b>HU</b>	Levenhuk MED 5M Mikroszkóp digitális kamerák 9,4" LCD képernyővel	<b>58</b>
<b>IT</b>	Levenhuk MED Videocamera digitale per microscopio da 5M con schermo LCD da 9,4"	<b>69</b>
<b>PL</b>	Levenhuk MED 5M Cyfrowa Kamera mikroskopowa z 9,4" LCD ekranem	<b>80</b>
<b>PT</b>	Levenhuk MED 5M microscópio digital câmera com ecrã LCD de 9,4"	<b>91</b>
<b>RU</b>	Цифровая камера Levenhuk MED 5 Мпикс с ЖК-экраном 9,4" для микроскопов	<b>102</b>


Never look at the sources of bright light or lasers through the camera or the microscope, as this may cause permanent eye damage and even blindness.

### General information

Levenhuk MED Digital Camera is a professional accessory for trinocular microscopes. Its 9.4" color LCD touch screen and 5Mpx digital camera allow you to conduct high-precision research in real time, and save images and video for further use.

### Parts of digital camera:

1. Microphone
2. USB port
3. USB mouse port
4. Headphone jack
5. AC adapter socket
6. MicroSD (TF card) slot
7. MiniHDMI out
8. Power on/off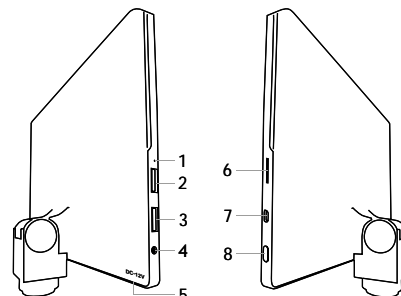

Caution! Remember that mains voltage in the U.S. and Canada is 110V and 220–240V in most European countries. Please refer to the specifications table for the correct mains voltage and never attempt to plug a 110V device into 220V outlet and vice versa without using a converter.

### Installing a digital camera

LCD screen digital camera is installed on the third vertical tube of trinocular microscopes. Assemble and set up your microscope. Unscrew the upper part of the trinocular tube and then insert (screw in) the camera. Connect the supplied AC adapter to the camera and to the AC network. Press the power button and the camera will be ready to work.

### Embedded software

The embedded application is designed for using Android digital cameras with microscopes. With this application, you can change image parameters, measure objects as well as save images and videos for further use. In addition, you can perform particle analysis with its image processing feature. Like all Android applications, this application features a user-friendly multilanguage interface.


## Camera display interface

The screen is divided into two parts:

1. The left part is the Preview Image. It displays the current image provided by the camera.
2. The right part is the Tool Panel. It contains a number of tools for image capturing, changing parameters, measuring, etc. Use the right side of the screen to perform the following actions:
  - Effect: Settings of the camera parameters
  - Measure: Measuring tool
  - Playback: Images and video playback
  - Snap: Image capturing
  - Record: Video recording
  - Setting: Setting and changing parameters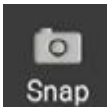

## Capturing and playback tools

 Snap	Press the Snap button to take a photo. Select the image size in the Settings section.
 Record	Press the Record button to start video recording. To stop recording, press the same button once again. If the original size of the video is smaller than 1080p, then the size of the recorded video will be the same as the original. If the original size is bigger than 1080p, the recorded video size will be reduced to a size smaller than 1080p.
 Playback	Press the <b>Playback</b> button to open the last image captured or to play the last video recorded. The image will be opened in the <b>Image Analysis</b> window.

## Camera settings

The **Effect** panel allows you to edit the captured image, if necessary.

### Exposure

Auto Exposure: If you enable the checkbox next to “Auto Exposure”, the brightness of the image will be automatically adjusted.

Manual mode: You can set the exposure time and gain.

### Color temperature (White balance)

Auto mode: If you enable the checkbox next to “Auto WB”, the white balance will be automatically set.

Manual mode: For manual settings, use the “Color temperature” slider tool.

**Brightness**: Use the slider tool to adjust the brightness.

**Contrast**: Use the slider tool to adjust the contrast.

**Saturation**: Use the slider tool to adjust the saturation.

**Sharpness**: Use the slider tool to adjust the sharpness.

**Gamma**: Use the slider tool for the gamma correction.

**Flip image**: These icons are used to flip the image vertically or horizontally.

This icon is for saving the settings or resetting all of the settings to the default parameters. You can save up to four different parameter sets as well as return to the factory settings.

## Measurements

On the Measure panel, you can select a tool to conduct a measurement.


Before you start, you need to calibrate the ruler. You also need to calibrate it after changing the microscope objective and/or camera resolution.

### 1. Calibrations

Go to the Measure panel, and press Add button to start a new calibration process. Press Edit button to make changes in previous calibrations.


Open calibration menu: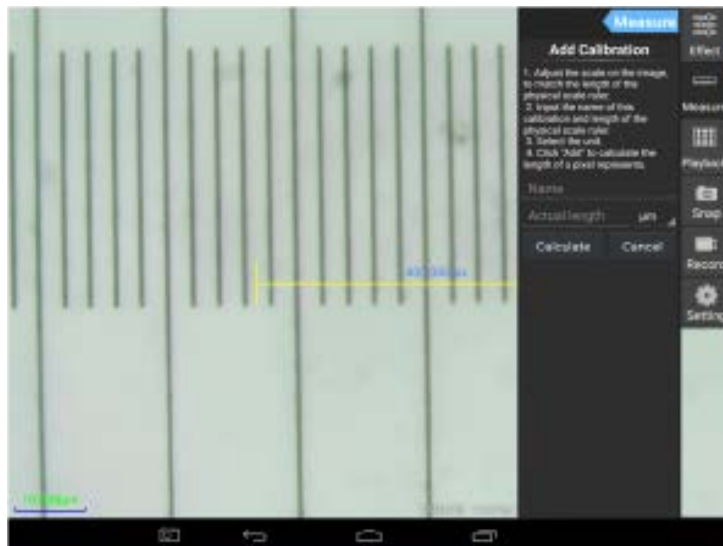


Follow the instructions:

Adjust the ruler to where its ends are on the physical ruler's scales. The ruler on the screenshot has 0.01mm scales, with each big grid equaling 10µm, and the whole ruler equals 1000µm.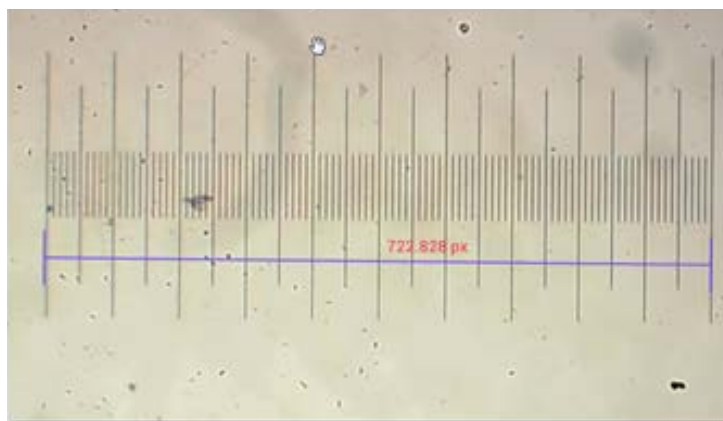

Enter the name for the calibration and actual length of the ruler. In the example, the name of the calibration is "4x", which corresponds to the magnification value of the chosen eyepiece, and 1000µm is the actual length of the chosen ruler.

Press Calculate to get the calibration value for the current eyepiece and image size. Press Add to save the calibration.

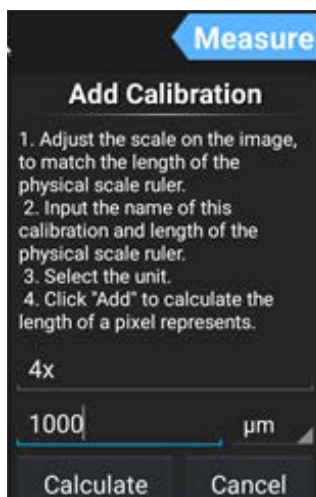
Calibrations		
Name	Data	Unit
4x	1.3842217	µm
10x	0.10938032	µm

Add Edit Delete
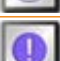
## 2. Measure tool

You can use a wide range of measure tools (rulers).

Icon	Measure tool	Action
	Point counting	Set the starting point of measuring on the photo
	Line	Measure the distance between two points
	Rectangle	Measure the width, length, and area of the rectangle
	Circle	Measure the area of a circle
	Cross	Measure the cross hair
	Angle	Measure the angle
	Two circles	Measure the distance between the centers of two circles
	Perpendicular	Measure the length of a perpendicular line
	Concentric circles	Measure the radius of two concentric circles
	Text annotation	Add a comment to an image
	Option	Change the width or the color of the ruler's scale, and the size or the color of the text
	Export	Export the image with measure marking
	Delete	Delete the chosen ruler
	Delete	Delete all


## 3. Use of the measuring ruler

Choose a calibration (the chosen option will be displayed as a black print on the white background).

Calibrations		
Name	Data	Unit
4x	1.3842217	µm
10x	0.10938032	µm


Add Edit Delete

Choose a ruler from the side menu, and it will appear in the image.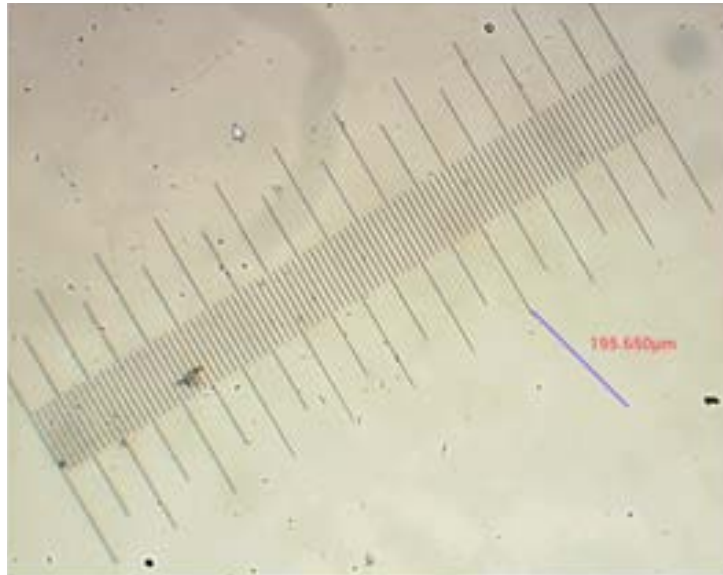

Use the measuring ruler to measure the physical ruler.

In this example, 6 large grids were measured: the measuring ruler shows 597.472µm, which confirms the correctness of the measurements.

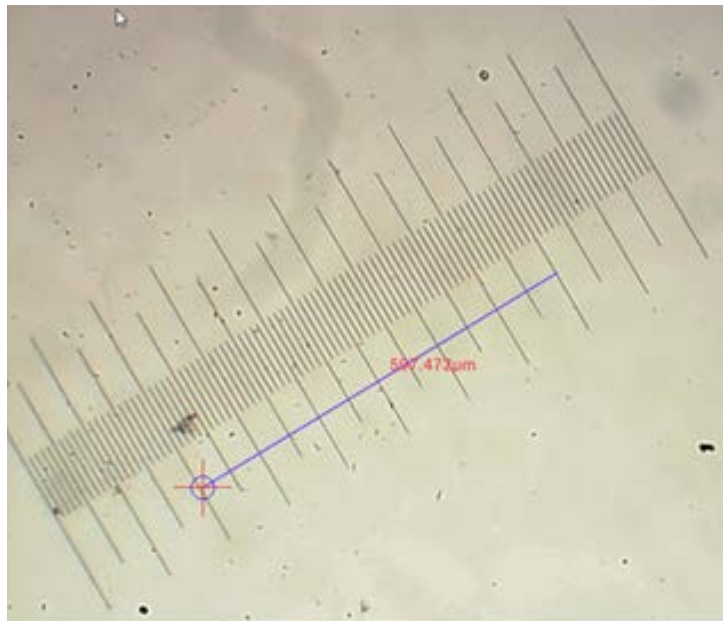
## Image analysis


The images are analyzed with the Android application YWImage, which uses OpenCV (Open Source Computer Vision Library). The application version 1.0 supports such functions as gray scale, contrast, binarization, and particle analysis.

To enable this feature, select the **Enable built-in image viewer** option in the settings dialog box.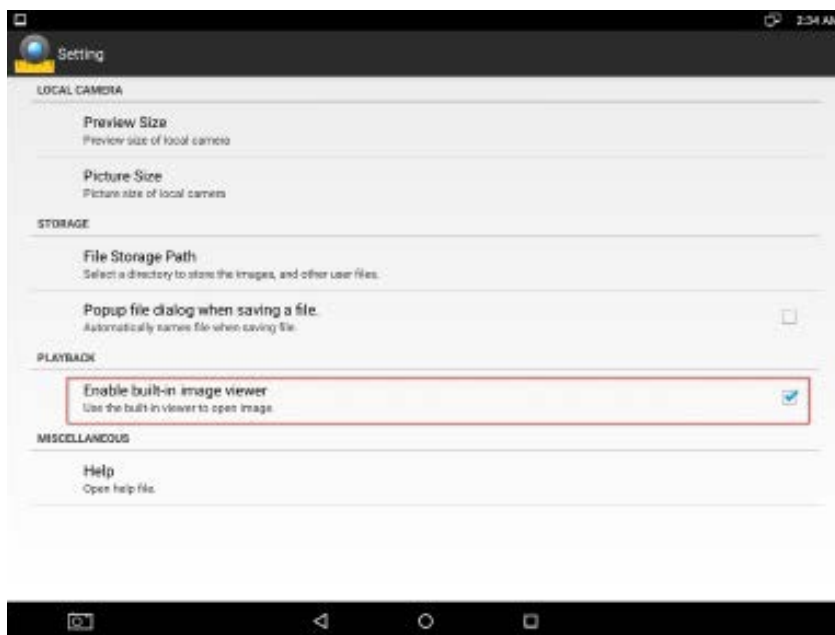


Let us consider an example of particle analysis.


Image uploading. Click the camera icon in the upper right corner of the screen. You can select an existing image or take a picture with the camera.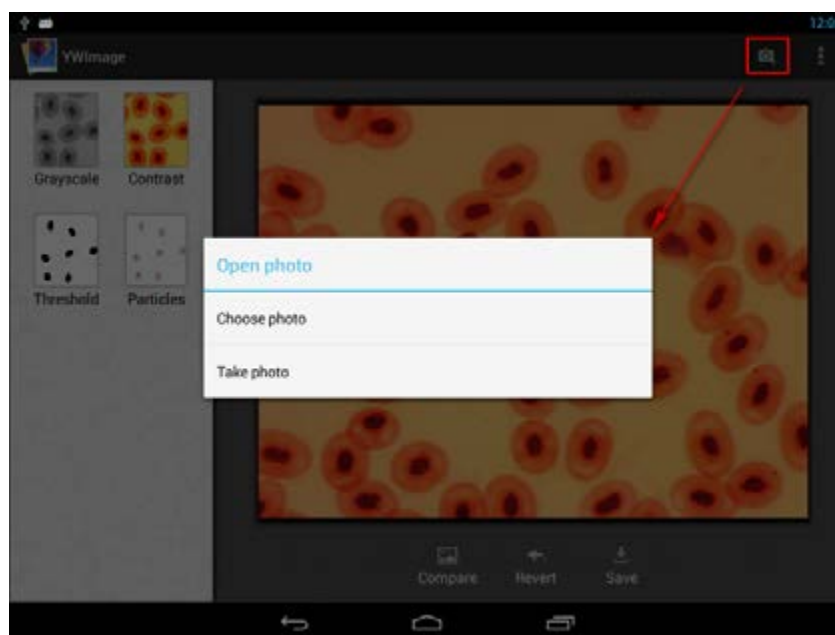

Convert a color photo to Grayscale mode. Choose **Grayscale** from the left menu and press **Apply** to confirm converting.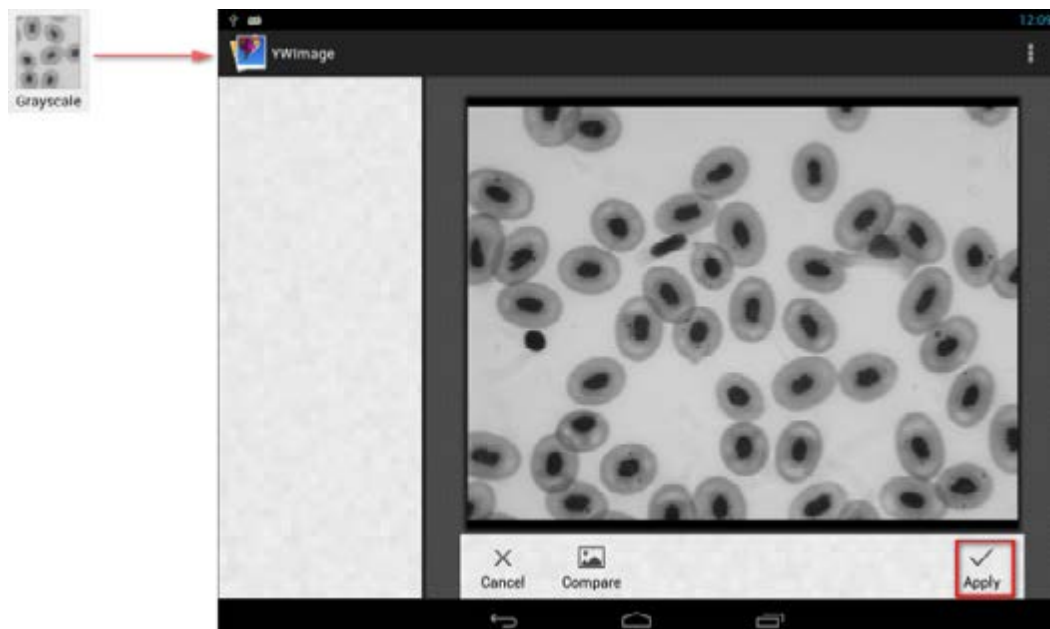


Contrast and brightness adjustment. Adjust the contrast and brightness of the black and white image, increasing the difference between objects and the background.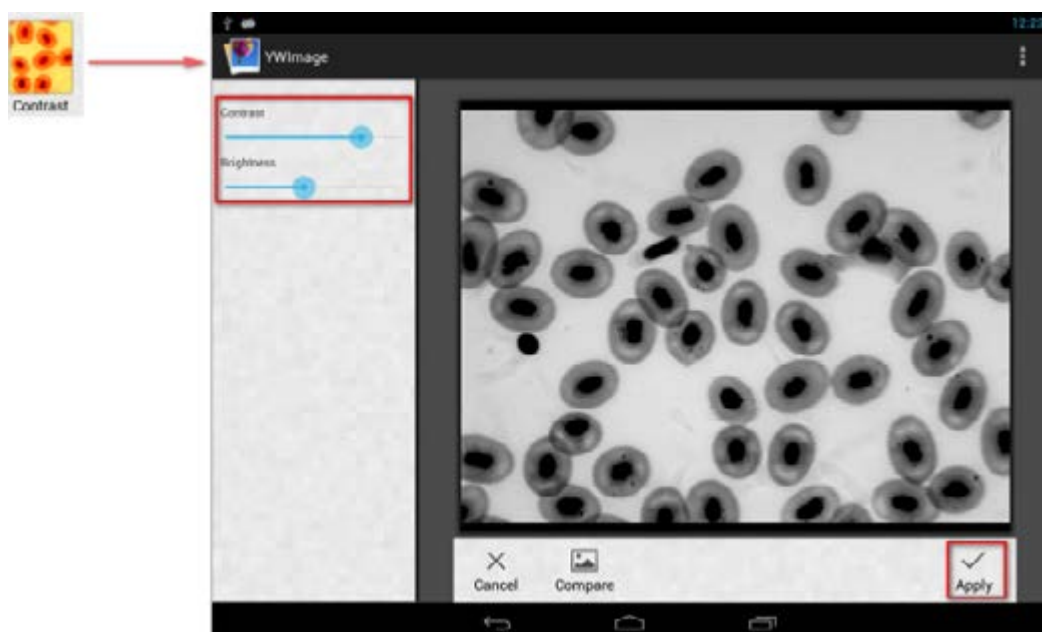

**Binarization.** Using the binarization function, you can convert a color image or an image in grayscale to two-color (black and white) by setting the minimum and maximum thresholds with which the brightness of each pixel will be compared.

All pixels within the threshold range will be highlighted in red. After executing the binarization, the RGB-value of the red areas will be set to (0, 0, 0), the RGB-value of the other sections will be (255, 255, 255).

**Note:** Red areas should not overlap.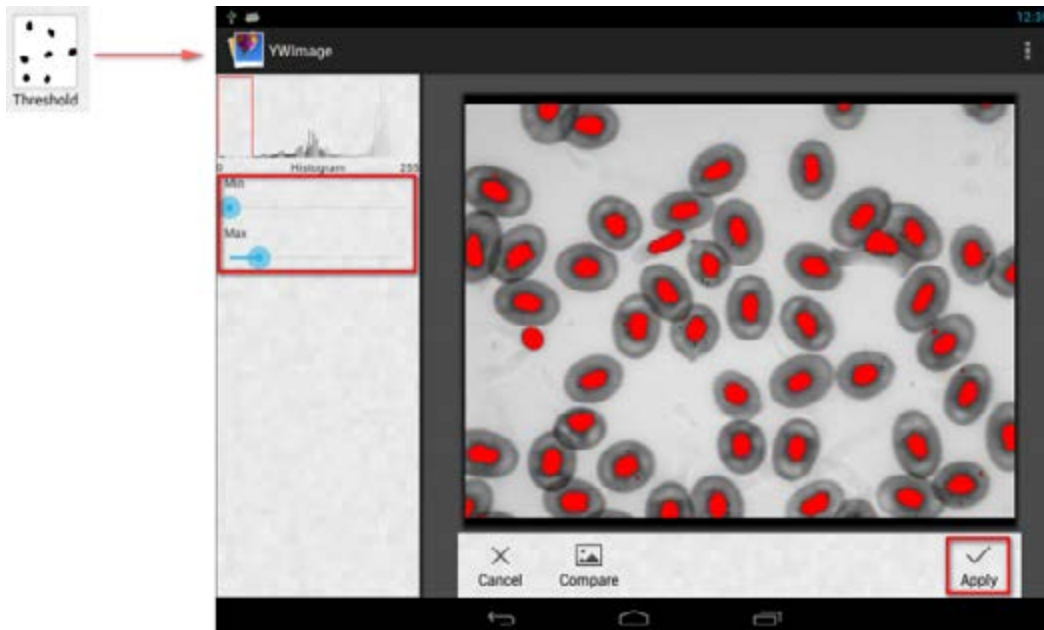


**Particle analysis.** After executing the binarization, open the particle analysis tool.

First of all, if some noises were not deleted during binarization, you need to specify a particle size range. The designation of the particle size range will make the result more accurate.

Press the button , to set the maximum value to infinity. Select the Show Outlines check box to show the boundaries and number of particles after the analysis. Click **Analyze Particles** to begin the analysis. The result will be shown below.

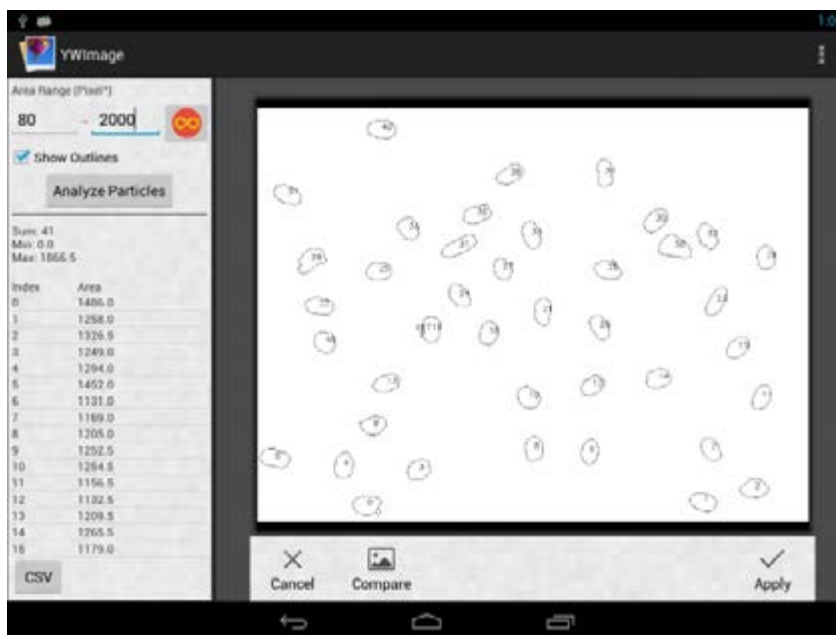Click the CSV button to export the data to the report file, which will be saved here by default: /mnt/sdcard/Image/Reports

## Specifications:

LCD display	9.4" color touch screen
Megapixels	5
Maximum resolution (for still images)	2048x1536px
Sensor	1/2.5"
Pixel size	2.2x2.2µm
Frame rate	15 frames per second
Exposure control	auto/manual
Video recording	+
Image format	*.jpg
Video format	1080p, *.3gp
White balance	auto/manual
Output	USB 2.0 (2pcs.), miniHDMI, Wi-Fi, TF card slot
Power supply	12V/2A; via AC adapter
Software features	measuring, brightness, exposure time, particle analysis, etc.
Mounting location	third eyepiece tube of a diameter of 23.2mm
OS	Android 5.1 (multilingual)
Embedded memory	4GB
Ability to connect additional equipment (not included in the kit)	support of microSD cards with capacity up to 32GB; monitor/TV (with a HDMI port); flash disk, mouse, keyboard (with a USB port); headphones (3.5mm)
The kit includes	digital camera with LCD screen, AC adapter

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Levenhuk reserves the right to modify or discontinue any product without prior notice.

## Care and maintenance

- **Never look at the sources of bright light or lasers through the camera or the microscope: it will cause PERMANENT DAMAGE TO YOUR EYES!**
- Do not disassemble the camera on your own.
- Keep the camera away from moisture; do not use it in rainy weather.
- Keep the camera away from shock or excessive pressure.
- Do not overtighten the locking screws.
- Keep the camera away from hazardous environment, home and car heaters, incandescent lamps or open fire.
- When cleaning the lenses, first blow any dust or debris off the surface or wipe them off with a soft brush. Then wipe the lens with a soft cleaning cloth slightly damp with spirits or ether.
- **Seek medical advice immediately if a small part or a battery is swallowed.**

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime warranty** against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk. This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc. For further details, please visit our web site: [www.levenhuk.com/warranty](http://www.levenhuk.com/warranty) If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk® is a registered trademark of Levenhuk, Inc.
© 2006–2021 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 928 E 124 Ave Ste D, Tampa, FL 33612 USA